

NEW! From Spectrum Laboratories, the Lab Dialysis Experts

Spectra/Por® Float-A-Lyzer® G2

Most Cost-Effective Ready-to-Use Dialysis Device

Spectrum introduces the next generation (G2) in Ready-to-Use laboratory dialysis devices featuring Ultra-pure Biotech Cellulose Ester (CE) Membrane.

- 95 – 98% Sample Recovery
- Highest Membrane Purity
- Superior Handling & Leak Protection
- Volume Specific Dilution Control
- Complete & Ready-to-Use Packaging
- Ideal for Gel Electro-elution

Sample Recovery, Purity and Dilution Control

Spectrum's proprietary Biotech Grade CE is a low protein-binding synthetic membrane available in 9 precise MWCO's with no heavy metal and sulfide contaminants. The cylindrical tubing geometry prevents sample dilution (associated with cassette-type devices) and provides open access for total volume retrieval. Only the Float-A-Lyzer G2 assures a 95-98% sample recovery while maintaining 99% sample purity and <1% sample dilution.

Float-A-Lyzer G2 - Cellulose Ester (CE) 12/pkg

MWCO	Color Code	Maximum Sample Volume		
		1 ml	5 ml	10 ml
0.1 - 0.5 kD	Green	G235025	G235049	G235061
0.5 - 1.0 kD	Orange	G235027	G235051	G235063
3.5 - 5 kD	Black	G235029	G235053	G235065
8 - 10 kD	Yellow	G235031	G235055	G235067
20 kD	Red	G235033	G235057	G235069
50 kD	Violet	G235034	G235058	G235070
100 kD	Blue	G235035	G235059	G235071
300 kD	Amber	G235036	G235060	G235072
1,000 kD	Pink	G235037	G235062	G235073



REQUEST FREE SAMPLE at SPECTRUMLABS.COM

Designed for Ease-of-Use and Leak Protection

The leak-proof screw-on cap with sealing o-ring provides easy access with included pipette for loading, in-process testing and sample retrieval, without the risk of needle punctures. The included flotation ring improves sample buoyancy and vertical orientation during dialysis. The sleek design allows multiple samples to be dialyzed in the same reservoir.

No additional kits, buoys or accessories are required. The Float-A-Lyzer G2 is packaged dry and ready-to-go with all components included.

- Simply rinse with water
- Load sample with Pipette
- And dialyze



go to website

NEW! Spectra/Por® Tube-A-Lyzer®

Disposable Ready-to-Use Dialysis Device for Rapid Separation and Dynamic Dialysis



Dialyze in 4-12 hours!

95 -98% Sample Recovery

2 Volume sizes (10 & 30 ml) & 6 MWCO's

Pyrogen-Free Biotech CE membrane for higher sample purity



go to www.spectrumlabs.eu for more info

